


ImageScope Upgrade:

- 1) Download the latest version of ImageScope: [Aperio ImageScope - Pathology Slide Viewing Software \(leicabiosystems.com\)](http://leicabiosystems.com)
- 2) Extract all files from the Zip folder
- 3) Run the .exe file and select “Remove”

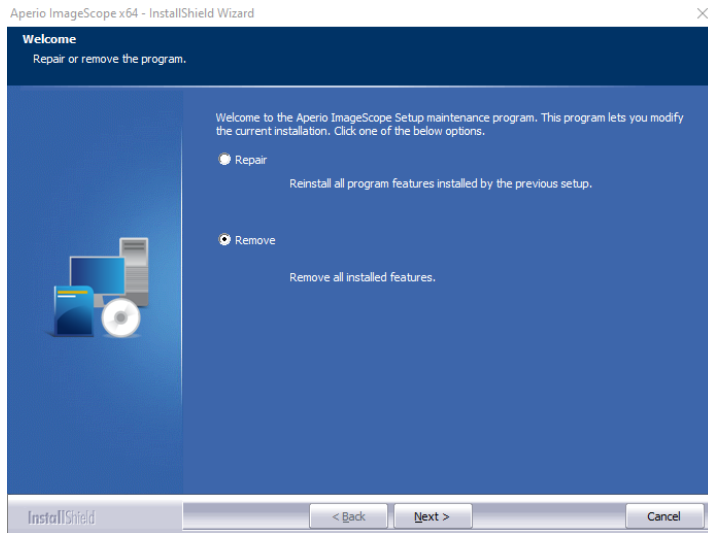

- 4) Restart your PC once finished
- 5) Go back to the .exe file and right click on it → Run as administrator (you may need assistance from your IT specialists)
- 6) Install program as recommended.

If ImageScope still doesn't work, do step #3 again but before restarting your PC, do the following steps:

- a) Open the Registry Editor (click on the Windows icon on the keyboard and type “Registry editor”)
- b) Go to: HKEY_CURRENT_USER → SOFTWARE → VB and VBA Program Settings → ImageScope64
- c) Right click on ImageScope64 → Delete
- d) Restart your PC
- e) Re-launch the .exe file
- f) Install the program as recommended

ImageScope should now work on your computer. If not, please reach out to Charley Assenmacher (chasse@upenn.edu). A “hotfix” version of ImageScope is available for the rare case where the downloaded version won't work.